



Chlorine-free formula tackles chemical spills  
in addition to standard blood / fluid spills.

# COMBAT SPILLS

*Solidify Chemical Spills, Blood, and  
Bodily Fluid Spills with Ease.*



## Green-Z® Spill Control Solidifier (Chlorine-Free Formula)

At times, workers may be unintentionally exposed to blood and bodily fluid spills along with other harmful substances such as toxic chemicals, biological hazards and chemotherapy agents. To help reduce those exposure risks, rely on Green-Z® for improved safety in your spill cleanup process.



**Fast-Acting**



**Chlorine-Free**



**No Splashing**

- **Contains fast-acting, superabsorbent polymers**
- **Chlorine-free** (safe to use on chemotherapy, glutaraldehyde, and mercury spills)
- **Turns spills into semi-solid (gel) mass**
- **Cost-effective cleanup solution**

**1oz** (approx. 30cc) : **1L** (1,000cc) \*



\*1oz solidifies 1L of water-based fluid

### Perfect for Unexpected Spills in:

- Healthcare & Laboratories
- Occupational Health & Safety
- Correctional Facilities
- Industrial & Manufacturing
- Transportation
- Schools
- Restaurants



**Safetec of America, Inc.**  
1-800-456-7077 | safetec.com

## Directions for Use

1. Wear appropriate protective clothing.
2. Sprinkle Green-Z powder to surround the spill and then evenly distribute over it. Allow sufficient time for solidification to occur.
3. Remove solidified material with scoop and scraper (or similar tool) and place in an appropriate container\*\*\*.
4. Clean and disinfect contaminated areas after use and wash thoroughly.



**Wear Protective Clothing**



**Sprinkle Green-Z® Over Spill**



**Remove Solidified Spill**

## Expiration Dating

Green-Z® is shelf stable and does not require an expiration date. In order for the product to remain shelf stable, it must be kept in a cool, dry place and in its original packaging. If contents are hard or compact, simply shake the product to loosen, before use.

\* Careful consideration should be given to chemical spill cleanup, i.e., room size, temperature, quantity and concentration(s) of the solution(s).

\*\*\* Dispose of infectious and hazardous waste in accordance with all local, state and federal regulations.

## Green-Z® Order Information




Order No.	Description	Case Qty.
42019	Green-Z® 21g single-use pouch	100/case
42010	Green-Z® 5 oz. shaker top bottle	24/case
42012	Green-Z® 15 oz. shaker top bottle	12/case

## Kits Containing Green-Z®



Order No.	Description	Case Qty.
49725	Chemotherapy Spill Kit (poly bag)	12/case
15201	Multi-Purpose Kit (hard case)	1 each
15205	Spill Leader Kit (hard case)	1 each

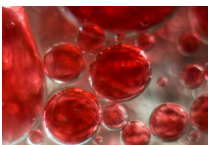

### OSHA Compliance 1910.1030 – Bloodborne Pathogens Standard

Universal Precautions is an approach to infection control. According to the concept of Universal Precautions, all human blood and certain human body fluids are treated as if known to be infectious.



**Safetec of America, Inc.**  
1-800-456-7077 | safetec.com